

---

---

## UNCP Emergency Eyewash/Safety Shower Weekly Inspection Checklist

**Prior to flushing the eyewash and safety shower stations, check for the following:**

- Is the area around the station clear of obstructions?
- Are the dust caps/covers over the eyewash nozzles in place?

**Please insure that a bucket is used to capture water from the station during the test and clean up any spilled water. Place a wet floor sign on the floor if required.**

**To test the eyewash:**

- Depress the paddle until water flows. The water should remain flowing after you remove your hand from the paddle.
- Continue to flow water until the water is clear with no indication of rust, dirt, etc. for 10-15 seconds.
- Feel the water to make sure it is tepid (room temperature - not hot or cold).

**To test the safety shower:**

- Pull the handle of the safety shower and flow water for 15-20 seconds. The water should flow without the handle having to be held.

**Room# \_\_\_\_\_ If any problems are found, e-mail the EH&S Office at [safety@uncp.edu](mailto:safety@uncp.edu).**

Year:	Week 1	Week 2	Week 3	Week 4	Week 5	Comments
January						
February						
March						
April						
May						
June						
July						
August						
September						
October						
November						
December						